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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5308-0350 / 5308-0700 / 5308-1050

AquaSampler

AquaSampler

AquaSampler

AquaSampler

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument is not suitable for sampling highly flammable/explosive liquids.

WARNING!

Damage to the respiratory tract and other organs due to escaping or outgassing media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Only use the instrument in safe environmental conditions.
- ▶ The sampling area has to be free of disturbing factors and hazard sources (e.g. rotating elements, currents, people).
- ▶ Secure the sampling area before sampling and ensure to have a secure footing.

- ▶ Do not store the instrument when it is filled.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The instrument is suitable to take samples of liquids and sediments.
- ▶ The instrument is suitable for sampling liquids from 0 - 60 °C (use at higher temperatures on request).
- ▶ The instrument is suitable to take samples up to 50 m depth (including working area).
- ▶ Minimum diameter of the sampling area: Diameter of the dipping bottle x 3.
- ▶ The instrument is only suitable to take samples and not for continuous operation or for storage.
- ▶ The instrument is maintenance-free.

Parts in contact with medium

- ▶ FEP/PTFE

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

Attach a suitable lowering cable (see accessories) securely to the eyelet ② at the top part of the Aqua-Sampler using a suitable cable clamp.

The AquaSampler can be extended using the extension tubes and the coupling (see accessories).

WARNING!

Bodily injury due to breakage or falling down during operation.

- ▶ Ensure correct handling.
- ▶ Ensure to attach the lowering hose securely to the immersion cylinder.
- ▶ Check all parts for a firm fit and damages before use.
- ▶ Wear the necessary protective clothing.

Safe sampling

1. Use a lowering cable with a suitable device (e.g. manually-operated reel (see accessories)) to lower the AquaSampler ① into the medium to be sampled. Do not lower the instrument in free fall and do not let it swing.
2. During lowering the liquids flows through the AquaSampler.

3. At the desired sampling depth pull the AquaSampler back using the lowering cable. Thereby, the bottom valve ③ closes automatically.

CAUTION!

Bodily injury due to the lowering cable slipping out of the reel or crushing of fingers.

- Pay attention to coil up lowering cable straight.
- Keep fingers away from cable guidance groove.

4. Pull the AquaSampler out of the medium to be sampled using the lowering cable.
5. Empty the sample through the spout ④ into a suitable collecting container.

WARNING!

Bodily injury due to escaping media when emptying the dipping bottle. Parts of the instrument can be filled with media.

- Empty the dipping bottle carefully.
- Wear the necessary protective clothing.

After taking the sample

Screw off the bottom valve and the spout and rinse all parts with a pH-neutral liquid.

Troubleshooting

Problem	To eliminate the error
Bottom valve is dripping.	Bottom valve is soiled. Screw off the bottom valve and clean it.

Accessories

Type	Specification	Item No.
Extension tube	350 ml	5308-0351
Extension tube	700 ml	5308-0701
Extension tube	1050 ml	5308-1051
Coupling	for extension tube	5308-0003
Manually-operated reel	PP	5309-0010
Stainless steel lowering cable, PTFE-coated	Ø 2 mm, Length 10 m	5309-0110
Lowering cable made of PP cord	Ø 2 mm, Length 10 m	5309-0201
Stainless steel lowering cable	Ø 1,25 mm, Length 10 m	5309-0301
The lowering cables are also available in lengths of 25 m and 50 m.		

More accessories can be found at www.buerkle.de

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