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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5330-1060 / 5330-1100 / 5330-1200 / 5330-2060 / 5330-2100 / 5330-2200

LiquiSampler

LiquiSampler

LiquiSampler

LiquiSampler

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

WARNING!

Damage to the respiratory tract and other organs due to escaping or outgassing media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

- ▶ The device can be used in the temperature range of 0 - 60 °C.
- ▶ Do not perform any lever movements during the application. The instrument may be deformed.

Layout and function

- ▶ The instrument is made of PP (item no. 5330-1xx0) or PTFE/FEP (item no. 5330-2xx0).
- ▶ The instrument is suitable to take samples of liquids.
- ▶ The instrument is only suitable to take samples and not for continuous operation.
- ▶ The instrument is not suitable for the storage of samples.
- ▶ The instrument is maintenance-free.

Parts in contact with medium

- ▶ PP (item no. 5330-1xx0)
- ▶ PTFE/FEP (item no. 5330-2xx0)

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Safe sampling

WARNING!

Bodily injury due to uncontrolled escaping liquid under pressure.

- ▶ Wear the appropriate protective clothing.

Point sampling/soil sample

1. Slide the wiping ring ① to the desired sampling depth.
2. Submerge the instrument into the medium to be sampled up to the desired depth / to the ground.
3. Push the push button ②. Thereby, the bottom valve opens and the instrument is filled with the liquid to be sampled.
4. Release the push button. Thereby, the bottom valve closes.
5. Remove the instrument from the medium to be sampled.
6. Empty the sample into a suitable collecting container by pushing the push button.

WARNING!

Bodily injury due to uncontrolled escaping liquid.

- ▶ Do not tilt the instrument when it is filled. Liquids may escape through the ventilation at the push button.
- ▶ Wear the appropriate protective clothing.

All-layer sample

1. Slide the wiping ring ① to the desired sampling depth.
2. Push the push button and hold it. Submerge the instrument into the medium to be sampled up to the desired depth. Thereby, the instrument is filled with the liquid to be sampled.
3. Release the push button. Thereby, the bottom valve closes.
4. Remove the instrument from the medium to be sampled.
5. Empty the sample into a suitable collecting container by pushing the push button.

After taking the sample

1. Slide the wiping ring with the aid of the supplied pusher down to wipe any remaining liquids.
2. Screw off the discharge funnel ③.
3. Hold the valve cone and screw off the push button.
4. Pull all parts out of the sampling tube.
5. Clean all parts. Use the cleaning brushes (see accessories) for the inside of tubes.

Troubleshooting

Problem	To eliminate the error
The instrument doesn't fit through the container opening.	Screw off the discharge funnel and take the sample without it. Screw the discharge funnel on again before emptying the sample.
Bottom valve is dripping.	<ol style="list-style-type: none"> 1. Bottom valve is soiled. Clean the instrument. 2. Discharge funnel is screwed on too tightly and the valve seat is deformed. Unscrew the discharge funnel.

Accessories

Type	Specification	Number	Item No.
Cleaning brush	PVC/stainless steel, 100 cm	1	5304-0100
Drain safety stand	Total length 117 cm, Ø 35 mm	1	5618-1000
Sample bottle HDPE	250 ml	1	0342-0250
Sample bottle HDPE	500 ml	1	0342-0500

More accessories can be found at www.buerkle.de

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