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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5365-6000

Tauchbombe Target

Target immersion cylinder

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Bombe à immersion Target

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument is not suitable for sampling highly flammable/explosive liquids.

WARNING!

Damage to the respiratory tract and other organs due to escaping or outgassing media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/disinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Only use the instrument in safe environmental conditions.
- ▶ The sampling area has to be free of disturbing factors and hazard sources (e.g. rotating elements, currents, people).
- ▶ Secure the sampling area before sampling and ensure to have a secure footing.
- ▶ Do not store the instrument when it is filled.

- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The instrument is suitable to take samples of liquids and sediments.
- ▶ The instrument is suitable for sampling liquids from 0 - 60 °C (use at higher temperatures on request).
- ▶ The instrument is suitable to take samples up to 5 m depth (including working area).
- ▶ Minimum diameter of the sampling area: Diameter of the dipping bottle x 3.
- ▶ The instrument is only suitable to take samples and not for continuous operation or for storage.
- ▶ The instrument is maintenance-free.

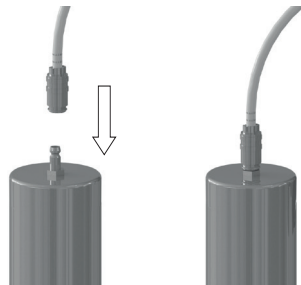
Parts in contact with medium

- ▶ Nickel-plated brass (immersion cylinder, rapid coupling)
- ▶ PE (lowering hose)
- ▶ PP (manually-operated reel)
- ▶ PTFE (sealing tape)
- ▶ Stainless steel V2A (hose clamp)

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

Connect the lowering hose ① to the immersion cylinder ③ using the rapid coupling ②.

WARNING!

Bodily injury due to breakage or falling down during operation.

- ▶ Ensure correct handling.
- ▶ Ensure to attach the lowering hose securely to the immersion cylinder.
- ▶ Check all parts for a firm fit and damages before use.
- ▶ Wear the necessary protective clothing.

Safe sampling

1. Close the valve ④.
2. Use the lowering hose and the manually-operated reel to lower the immersion cylinder into the medium to be sampled. Do not lower the instrument in free fall and do not let it swing.

CAUTION!

Bodily injury due to the lowering hose slipping out of the reel or crushing of fingers.

- Pay attention to coil up lowering hose straight.
- Keep fingers away from cable guidance groove.

3. At the desired sampling depth open the valve. Thereby, the immersion cylinder is vented and filled.
4. Close the valve.
5. Pull the immersion cylinder out of the medium to be sampled using the lowering hose.
6. Disconnect the lowering hose from the immersion cylinder using the rapid coupling.
7. Screw off the top part ⑤ of the immersion cylinder.
8. Empty the sample into a suitable collecting container.

WARNING!

Bodily injury due to escaping media when emptying the dipping bottle. Parts of the instrument can be filled with media.

- Empty the dipping bottle carefully.
- Wear the necessary protective clothing.

After taking the sample

Screw off the top part and rinse the immersion cylinder and the lowering hose with a pH-neutral liquid.

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