## Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

## Safety instructions

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.

### WARNING!

When using with flammable liquids, do not spray into ignition sources.

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The device is suitable for indoor and outdoor use.

### CAUTION!

Consider possible influences of wind or draught during spraying.

- ▶ Wear the necessary protective clothing.

- ▶ The instrument is not suitable for cleaning in a dishwasher.

## Layout and function

- ▶ The device is made of PE, PP, FKM and V4A.
- ▶ The device is suitable for spraying particle-free liquids. The device is not suitable for continuous operation.
- ▶ The device can be used for liquids from 0 - 30 °C and can be used in the temperature range from 5 - 35 °C. Avoid direct sunlight.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

## Safe usage

1. Fill the bottle ① with liquid up to the „MAX“ mark.

### CAUTION!

An even and perfect atomization depends on the medium filled.

- ▶ Mix media thoroughly before filling.
- ▶ Do not overfill the container.

2. Screw spray head ② and tighten hand-tight.
3. Use pump rod ③ to pump air into the bottle (approx. 20 to 30 pumping movements).

### CAUTION!

The operating pressure is max. 4 bar.

- ▶ If the pressure is too high, the built-in safety valve opens automatically.

4. Press spray trigger ④ for spraying.

### WARNING!

Bodily injury due to uncontrolled spraying of fluid.

- ▶ Wear the necessary protective clothing.

5. Turn the nozzle nut ⑤ to adjust the spray jet. (From fine mist to powerful jet.)

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1. Release spray trigger ④ after spraying.
2. Open the spray head slowly to let escape the pressure.

### CAUTION!

Always let escape the pressure.

- ▶ Do not leave the device under pressure.

3. Clean the device.

### CAUTION!

Oil the pump head from time to time (see arrow on the spray button).

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