Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

! WARNING!

The instrument is not suitable for highly flammable or explosive liquids.

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.


Safety instructions during and after use

- ▶ The instrument generates overpressure. Possible bodily injury due to escaping liquids.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The instrument consists of a riser pipe made of polypropylene (PP), a discharge tube made of PVC, a shut-off valve made of PP and PE and a universal bung made of PVC.

- ▶ The instrument is suitable to fill off highly fluid and low-viscosity liquids.
- ▶ The bung fits onto container openings with a diameter of 30 to 70 mm forming an airtight seal.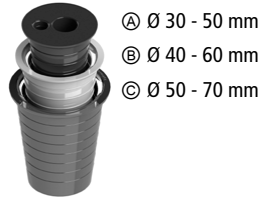

- ▶ The instrument is maintenance-free.

Point out any damage in the packaging to the freight forwarder/supplier immediately.


Assembly

! WARNING!

Some liquid will continue to run while the discharge hose is being screwed on. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve on the discharge hose before mounting.


1. Pull out the telescopic riser pipe ① completely.
2. Guide the riser pipe ① with the universal bung ② into the container to form an airtight seal.
3. If necessary adjust the universal bung to the opening diameter of the container.
4. Close the shut-off valve ⑥ on the discharge hose.
5. Plug the hand pump ball ④ into the bung.
6. Close the stop-valve ⑤ (push the lever upwards).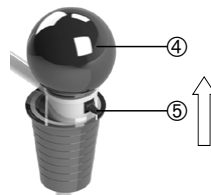

Safe filling

1. Hold the discharge hose in the filling container.
2. Activate the hand pump ball and generate overpressure. Liquid starts to flow.
3. Open the shut-off valve on the discharge hose.

4. To complete the filling process open the stop-valve (push the lever downwards) and allow the overpressure to escape or close the shut-off valve on the discharge tube.

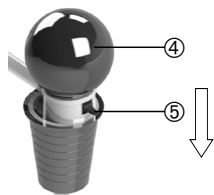**NOTE!**

The stop-valve is not suitable for precise dosing.

5. After filling is complete, close shut-off valve on the discharge hose.
6. To prevent any liquid continue to run (siphoning principle), fix the discharge hose above the liquid level in the original container.

Completing filling and dismantling**CAUTION!**

The container is under pressure. Bodily injury due to escaping liquid.

- ▶ Slowly remove the instrument from the container.

! WARNING!

Some liquid will continue to run during the removal of the discharge hose. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve on the discharge hose.
- ▶ Raise the discharge hose and allow the liquid to run back into the container.

1. Remove the hand pump ball from the bung.
2. Carefully remove the instrument from the container.
3. Rinse the instrument with a pH-neutral fluid.

Troubleshooting

| Problem | To eliminate the error |
|--------------------------------------------------|--------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------------|
| The instrument does not generate an overpressure | <ol style="list-style-type: none"> 1. Stop-valve is opened. 2. The instrument is not connected with the container to form a seal. |
| Instrument is not pumping any liquid | <ol style="list-style-type: none"> 1. The instrument is not connected with the container to form a seal. 2. The riser pipe does not extend into the liquid being filled. |

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