Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.
- ▶ The instrument is designed for use in industry, laboratories and science. The instrument is not suitable for any application in private or public spaces.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in well ventilated rooms or under a suitable exhaust hood. Wear appropriate protective clothing.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

WARNING!

Possible sparking due to electrostatic discharge. Danger of explosion when using combustible liquids.

- ▶ Connect the earthing terminal before use (see accessories).
- ▶ Wear the necessary protective clothing.

Safety instructions during and after use

- ▶ The instrument is suitable at a temperature range from 0 to 60 °C.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The instrument is made of stainless steel V2A, the seal is made of ePTFE.
- ▶ The instrument is to attach Bürkle barrel connectors and manual pumps (immersion tube Ø 32 mm) in vertical position in barrels with connection thread G 2".
- ▶ For container with nozzle (height of the container opening: 6-55 mm) and suitable for container openings of 48-75 mm diameter.

- ▶ The instrument is suitable for the gas-tight filling of liquids and slightly viscous media in conjunction with a manual pump (see accessories).

WARNING!

Possible escaping of liquid and gas due to excess pressure in the barrel. Bodily injury due to escaping gas and liquid.

- ▶ Avoid excessive excess pressure. (Do not expose the barrel to direct solar irradiation or variations in temperature. Pay attention to chemical resistance.)
- ▶ Ensure adequate air ventilation.
- ▶ Wear the necessary protective clothing.

NOTE!

Use a suitable pump to ensure gas-tight filling.

- ▶ It is recommended to use a pump that can be hermetically sealed by shut-off valves (see accessoires).

Scope of delivery

1 x	Universal adapter incl. ePTFE seal	①
1 x	Mounting key	②

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

1. With the mounting key ② screw the clamps ③ at the clamping jaws ④ in alternatel downwards (anti-clockwise) and turn them inward.
2. Place the thread adapter ① on the drum opening.
3. With the mounting key ②, screw the clamps ③ at the clamping jaws ④ alternately upwards (clockwise). Clamps ③ automatically turn outwards. Fix them hand-tight.

NOTE!

Clamps ③ must not block the opening.

- ▶ Point the clamps ③ outwards.

4. Mount the required pump or barrel connector.

NOTE!

Always hold the mounting key ② during assembly/dismounting to avoid possible loss of the mounting key ② in the container.

NOTE!

Ensure to use a fitting immersion tube (immersion tube Ø 32 mm, see accessories). Fix the immersion tube sufficiently.

NOTE!

Check correct connection.

- ▶ Screw on the universal thread adapter ① only hand tight and without force.

NOTE!

Regularly check seal for correct connection and integrity.

Dismounting

1. Disassemble the assembled pump or barrel connector.
2. With the mounting key ②, screw the clamps ③ at the clamping jaws ④ alternately downwards (anti-clockwise) and turn them inward.
3. Carefully remove the thread adapter ① from the container opening.
4. Rinse the thread adapter ① with a pH-neutral fluid.

CAUTION!

During disassembly of the thread adapter ①, medium may stick to the surface.

- ▶ Disassemble the pump and thread adapter ① carefully.
- ▶ Wear the necessary protective clothing.

Accessories

Type	Specification	Number	Item No.
PP barrel screw joint	R 2", steel fine outer thread	1	5600-3130
Brass barrel screw joint	R 2", steel fine outer thread	1	5601-0100
Gastight barrel connector	R 2", steel fine outer thread	1	5601-0130
PP barrel pump	Discharge hose/stopcock, Immersion depth 80 cm	1	5600-0802
Stainless steel barrel pump	Discharge tube, Immersion depth 91 cm	1	5601-0800
Stainless steel barrel pump	Discharge hose/stopcock, Immersion depth 91 cm	1	5601-0801
Solvent pump foot operated	Discharge hose/stopcock, Immersion depth 95 cm	1	5603-3000
Solvent pump hand operated	Immersion depth 60 cm	1	5603-1000
Withdrawal system for solvents	Discharge tube, Immersion depth 95 cm	1	5603-5003
Withdrawal system for solvents	Discharge hose, Immersion depth 95 cm	1	5603-5004
PTFE barrel pump Ultrapure	Discharge tube, Immersion depth 95 cm	1	5606-1000
PTFE barrel pump Ultrapure	Discharge hose, Immersion depth 95 cm	1	5606-1001
Gas-tight stainless steel barrel pump	Discharge hose, Immersion depth 91 cm	1	5611-0090
Anti-static set	To prevent electrostatic charging	1	5602-1000

More accessories can be found at www.buerkle.de

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