Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.
- ▶ The instrument is designed for use in industry, laboratories and science. The instrument is not suitable for any application in private or public spaces.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to escaping or off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood. Wear appropriate protective clothing.
- ▶ When working with explosive substances, follow the technical rules for combustible liquids (in accordance with TRBF – (German) Technical Rules on Combustible Liquids) and wear appropriate protective equipment (zone 0 and 1, see annex).

WARNING!

Possible sparking due to electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing terminal before use (see accessories).
- ▶ Wear the necessary protective clothing.

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Not suitable for acetone.
- ▶ Do not use any force when operating the instrument.

CAUTION!

Escaping liquids due to falling over or tipping of the instrument. Bodily injury due to escaping liquid and damage to the instrument.

- ▶ Ensure sufficient stability of the instrument.
- ▶ Shut down and secure the instrument immediately.
- ▶ Contact the dealer or manufacturer.

Safety instructions during and after use

- ▶ The instrument is suitable for indoor use and manual operation.
- ▶ Pump with manual operation. Not suitable for continuous operation.
- ▶ Pay attention to have sufficient space above during the filling process (minimum height: height of the maximum extended piston + 20 cm).
- ▶ The instrument is not suitable for cleaning in a dishwasher.

NOTE!

Pump and valves can be damaged or plugged by dirt particles.

- ▶ Do not fill off any polluted media.
- ▶ Rinse the pump with a pH-neutral medium.

Layout and function

- ▶ The instrument is made of conductive material (stainless steel 1.4301).
- ▶ The instrument is suitable for filling combustible liquids in accordance with groups IIA and IIB (see annex).
- ▶ The instrument is suitable for filling from containers with a capacity of 60 - 1000 l into smaller containers.
- ▶ The instrument is maintenance-free and suitable at a temperature range from 0 to 40 °C.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

WARNING!

Possible electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing terminal before use (see accessories).
- ▶ Wear the necessary protective clothing.

Assembly

WARNING!

Some liquid will continue to run while the discharge hose is being screwed on. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve on the discharge hose before mounting.

1. Screw the discharge hose ② onto the riser pipe ①.
2. Guide the barrel screw connection (accessory; not included in delivery) and riser pipe ① with the discharge hose ② into the container and screw it together finger-tight.
3. If necessary adjust the riser pipe.

NOTE!

Ensure to use a fitting thread (see accessories).

- ▶ Screw the thread by hand and without any force.

Safe filling

1. Hold the discharge hose in the filling container.
2. Open the shut-off valve ③ on the discharge hose.
3. Activate the piston rod ④ by using the handle ⑤ and fill off the liquid.

CAUTION!

A too heavy lateral load can deform the piston rod and cause a leakage of the pump head.

- ▶ Move the piston rod up and down in a straight line.

CAUTION!

Fluid ejection off the filling container. Bodily injury due to escaping liquid.

- ▶ Wear the necessary protective clothing.

Completing filling and dismounting

1. After the last pump stroke, leave the piston rod in the lower position and insert it into the pump up to the stop.
2. Close the shut-off valve on the discharge hose.
3. To prevent any liquid continue to run (siphoning principle), fix the discharge hose above the liquid level in the original container.

! WARNING!

Some liquid will continue to run after the discharge hose is unscrewed. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve on the discharge hose.
- ▶ Raise the discharge hose and allow the liquid to run back into the container.

4. Unscrew the discharge hose.
5. Carefully unscrew the instrument from the container.
6. Rinse the instrument with a pH-neutral fluid.

Accessories

Type	Specification	Number	Item No.
Antistatic set	To prevent electrostatic charging	1	5602-1000
Brass barrel screw joint	R2" steel fine outer thread	1	5601-0100
Gastight barrel connector	R2" steel fine outer thread	1	5601-0130
Discharge tube	For stainless steel barrel pump	1	5601-0080
Drain safety stand	Total length 117 cm, Ø 35 mm	1	5618-1000

More accessories can be found at www.buerkle.de

Troubleshooting

Problem	To eliminate the error
Liquid is not pumped.	Bottom valve is polluted. Clean the pump. If there is only little residual liquid in the container, the pumping must be carried out for a longer time.

Annex

Operational directive in accordance with 1999/92/EC:

Zone 0	a potentially explosive atmosphere is assumed to be continuously present
Zone 1	a potentially explosive atmosphere is assumed to be occasionally present
Zone 2	a potentially explosive atmosphere is not anticipated (and if it does occur, then only rarely and briefly)

Product directive in accordance with 94/9/EC:

Group I	for use in mining
Group II	for use in other areas subject to explosion hazard
Group IIA	Maximum permitted gap > 0.9 mm
Group IIB	0.5 mm < maximum permitted gap < 0.9 mm
Group IIC	Maximum permitted gap < 0.5 mm

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