**Bürkle GmbH**

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

## Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood.
- ▶ When working with explosive substances, follow the technical rules for combustible liquids (in accordance with TRBF – (German) Technical Rules on Combustible Liquids) and wear appropriate protective equipment (zone 0, see annex).

### ⚠ WARNING!

Possible sparking due to electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing terminal before use (see accessories).
  - ▶ Wear the necessary protective clothing.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

### NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

## Safety instructions during and after use

- ▶ The instrument generates overpressure. Possible bodily injury due to escaping liquids. Hold the shut-off valve on the discharge hose closed.
- ▶ The instrument generates overpressure. If there are differences in pressure, loud popping noises can be heard with some types of containers.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

## Layout and function

- ▶ The instrument is made of conductive material (1.4301).
- ▶ The instrument is suitable for filling combustible liquids in accordance with groups IIA and IIB (see annex).
- ▶ Always use the instrument with pressure-resistant containers (up to 220 l, >1 bar) for filling into open containers (<1 l).
- ▶ When filling combustible liquids, always use an antistatic set (see accessories).
- ▶ The instrument is maintenance-free and suitable at a temperature range from 0 to 60 °C.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

### ⚠ WARNING!

Possible electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Connect the earthing terminal before use (see accessories).
- ▶ Wear the necessary protective clothing.

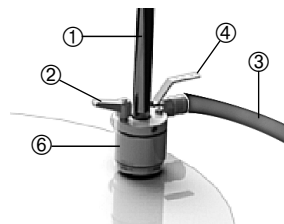
### ⚠ WARNING!

Some liquid will continue to run while the discharge hose is being screwed on. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve on the discharge hose before mounting.

## Assembly

1. Guide the barrel screw connection ⑥ and riser pipe ① into the container and screw it together finger-tight.
2. If necessary open the clamping lever ② and adjust the riser pipe. Then close the clamping lever and lock the riser pipe in place.
3. Screw the discharge hose onto the riser pipe ①
4. Attach the connecting hose to the foot bellows on the instrument ③.

### Safe filling

1. Open the shut-off valve ④ on the instrument.
2. Activate the foot bellows and generate overpressure.
3. Hold the discharge hose in the filling container.
4. Open the shut-off valve ⑤ on the discharge hose.
5. After filling is complete, close shut-off valve on the discharge hose.

#### **WARNING!**

The container is under pressure. Bodily injury due to escaping liquid.

- ▶ Move the shut-off valve on the instrument to a 45° position ④.
- ▶ Allow excess air to escape completely from the container.

#### **WARNING!**

Some liquid will continue to run after the discharge hose is unscrewed. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve on the discharge hose.
- ▶ Raise the discharge hose and allow the liquid to run back into the container.

### Completing filling and dismantling

1. Close the shut-off valve ④ on the instrument.
2. Unscrew the discharge hose.
3. Unscrew the connecting hose to the foot bellows ③.
4. Carefully unscrew the instrument from the container.
5. Rinse the instrument with a pH-neutral fluid.

### Accessories

Type	Specification	Number	Item No.
Antistatic set	to prevent electrostatic charging	1	5602-1000

More accessories can be found at [www.buerkle.de](http://www.buerkle.de)

### Annex

Operational directive in accordance with 1999/92/EC:

Zone 0	a potentially explosive atmosphere is assumed to be continuously present
Zone 1	a potentially explosive atmosphere is assumed to be occasionally present
Zone 2	a potentially explosive atmosphere is not anticipated (and if it does occur, then only rarely and briefly)

Explosion groups:

Group I	for use in mining
Group II	for use in other areas subject to explosion hazard
Group IIA	Maximum permitted gap > 0.9 mm
Group IIB	0.5 mm < maximum permitted gap < 0.9 mm
Group IIC	Maximum permitted gap < 0.5 mm

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