Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood.
- ▶ When working with explosive substances, follow the technical rules for combustible liquids (in accordance with TRBF – (German) Technical Rules on Combustible Liquids) and wear appropriate protective equipment (zone 0, see annex).

WARNING!

Possible sparking due to electrostatic discharge. Danger of explosion when using combustible liquids.

- ▶ Before use, insert the earthing cable connector ④ and connect the earthing terminal ②.
 - ▶ Wear the necessary protective clothing.
- ▶ Remove hazard sources within a radius of 50 cm around the instrument.
 - ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

Safety instructions during and after use

- ▶ Liquid may escape while the instrument is being unscrewed. Danger of bodily injury due to escaping liquid.
- ▶ The instrument must only be transported unfilled and without a container. Danger of bodily injury due to escaping liquid.

Layout and function

The instrument is made of conductive material (1.4301/1.4404). The instrument works according to the overpressure principle and is suitable for filling combustible liquids in accordance with group IIC (see annex). Always use the instrument with pressure-resistant containers (1 to 5 l, >1 bar) for filling into open containers (<1 l). The instrument is not suitable for cleaning in a dishwasher. The instrument is maintenance-free and can be used in the temperature range from 5 to 40° C.

Scope of delivery

| | | |
|-----|--|---|
| 1 x | Mini solvent pump | ① |
| 1 x | Earthing cable, length 2 m | ② |
| 1 x | PTFE hose (internal x external diameter: 6 x 8 mm), length 0.4 m | ③ |

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Mounting and safe filling

WARNING!

Possible electrostatic discharge. Danger of explosion if combustible liquids are used.

- ▶ Before use, insert the earthing cable connector ④ and connect the earthing terminal ②.
- ▶ Wear the necessary protective clothing.

CAUTION!

The metal thread of the instrument can damage glass containers. Bodily injuries due to glass splinters and escaping liquid.

- ▶ Hold the instrument vertical and screw it on carefully.

1. Attach the PTFE hose securely to the suction nozzle of the instrument ③. The PTFE hose must extend into the liquid being filled (shorten the PTFE hose if necessary and adapt it to the container you are using).
2. Screw the instrument onto the container air-tight.
3. Activate the hand pump ball manually.
4. Hold the filling container under the discharge and press the discharge button ⑤. If necessary a PTFE hose can be connected to the discharge (max. length 0.4 m).

CAUTION!

The container is under pressure. Bodily injury due to escaping liquid.

- ▶ Slowly unscrew the instrument from the container and allow the excess pressure to escape.

After filling

1. Carefully unscrew the instrument from the container.
2. Remove the PTFE hose.
3. Rinse the PTFE hose instrument with a pH-neutral liquid.

Troubleshooting

| Problem | To eliminate the error |
|--|--|
| The instrument does not generate an overpressure | 1. The instrument is not connected with the container to form a seal. |
| Instrument is not pumping any liquid | 1. The instrument is not connected with the container to form a seal. 2. The uptake hose is not connected with the instrument to form a seal. 3. The uptake hose does not extend into the liquid being filled. |

Annex

Operational directive in accordance with 1999/92/EC:

| | |
|--------|--|
| Zone 0 | a potentially explosive atmosphere is assumed to be continuously present |
| Zone 1 | a potentially explosive atmosphere is assumed to be occasionally present |
| Zone 2 | a potentially explosive atmosphere is not anticipated (and if it does occur, then only rarely and briefly) |

Explosion groups:

| | |
|-----------|--|
| Group I | for use in mining |
| Group II | for use in other areas subject to explosion hazard |
| Group IIA | Maximum permitted gap > 0.9 mm |
| Group IIB | 0.5 mm < maximum permitted gap < 0.9 mm |
| Group IIC | Maximum permitted gap < 0.5 mm |

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